

EyeSpray eye/face wash

Stainless steel

Article 850 1003

List of supplies

- EyeSpray rack stainless steel.
 - 2 spray bottles containing a sterile buffered sodium chloride solution (0.9 %).
 - Green and yellow instruction labels.
 - Red sealing tape and a white label for indication of expiry date.
 - Installation instructions (hereafter).
 - Instructions for replacement of the spray bottles (to be kept safe).
-

Installation instructions

- Fix carefully the green and yellow labels on the rack:
 - User instructions on the front of the rack
 - Bottle replacement instructions inside the rack
 - Above the handle (yellow label).
 - Determine a visible spot, close to the danger area.
 - Install the rack on the wall (screws not supplied), using if necessary the distance plastic washers if the wall is not flat.
 - The rack should be installed so that the spray bottles are at the height of 1.1 m, as for any eyewash, which corresponds to a distance of 1.15 m between the bottom of the closed rack and the floor.
 - Install the bottles according to enclosed replacement instructions.
 - Close the cover of the rack, leaving the red flag inside.
 - Fix the red sealing across the cover and the baseplate in such a manner that it will necessarily be broken when activating the eyewash. Also fix the white label with the expiry date.
 - These instructions can be found and a video can be seen on www.eyespray.ch
-

For questions, an order for new racks, replacement bottles or accessories, please contact your vendor, your plumber or directly the importer:

LaboSafe SA
Chemin du Petit-Pontarlier 15
CH – 2000 Neuchatel

☎ +41 32 756 96 96

www.labosafe.ch

info@labosafe.ch

EyeSpray eye/face wash

Stainless steel

Articles 850 1003

Replacement of the spray bottles

After each use, new bottles must be installed and the seal must be replaced. The red flag indicates that the eyewash has been activated and that new bottles must be installed.

- Open the cover of the rack, however not completely, leaving a space using a pen or an object thus allowing the spray bottles not to be activated.
- Dismount the bottles from their casing and remove them from the rack.
- Make sure to have two new bottles available, as well as a white label indicating the expiring date and a red sealing sticker.
- Write down on the white label the expiry date as indicated at the bottom of the bottle. The bottles are to be replaced after three years.
- Remove the cap from the new bottles.
- Install the new bottles in the rack, making sure the spray is facing upwards and the bottles are firmly fixed in their casing. A sticker on the cap of the bottle reminds about the need to correctly orient the nozzles.
- Close the cover of the rack, leaving the red flag inside, thus showing the rack has not been opened after replacement of the bottles.
- Fix the red sealing across the cover and the baseplate in such a manner that it will necessarily be broken when activating the eyewash. Also fix the white label with the expiry date, preferably on the side of the rack.

A periodical control is necessary to make sure that the seal is not broken, that the red flag is not visible and that the expiry date has not been reached.

These instructions can be found and a video can be seen on www.eyespray.ch

For questions, an order for new racks, replacement bottles or accessories, please contact your vendor, your plumber or directly the importer:

LaboSafe SA
Chemin du Petit-Pontarlier 15
CH – 2000 Neuchâtel

☎ +41 32 756 96 96

www.labosafe.ch

info@labosafe.ch